



S-Pak® Membranes

An easy-to-use membrane filter

To ensure maximum productivity and efficiency, each individually packed S-Pak® membrane is protected with a blue sheet of paper that keeps the membrane from coming into contact with the film. This protective sheet prevents the migration of any adhesive.

S-Pak® Membrane Advantages

- Individually packed
- Sterile
- · Ready to use
- Easy to use
- No hydrophobic spots
- No toxicity
- Better recoveries





Easy to use and manipulate with gloved hands



Utilizing sterile forceps, peel back the film and separate the blue paper from the membrane. Remove the membrane with the forceps.



Place the membrane and blue paper on a wet filtration support. When wet, the blue paper curls enabling you to easily remove it using your forceps.

Ordering Information

S-Pak® Membrane Filters: 600 (4 boxes of 150 membranes)

Cat. No.	Filter Color	Filter Diameter	Dava Ciza (um)
Cat. No.	Filter Color	Filter Diameter	Pore Size (µm)
GSWG047S6	white	47mm	0.22
HABG047S6	black	47mm	0.45
HAWG047S6	white	47mm	0.45
HVWP047S6	white	47mm	0.45 (PVDF)
HCWG047S6	white	47mm	0.7
AABG047S6	black	47mm	0.8
AAWG047S6	white	47mm	0.8
RAWG047S6	white	47mm	1.2

We offer a range of membranes to meet your specific filtration requirements. Please contact your local sales representative for additional information.

For your high volume applications, try our EZ-Pak® membrane filters!

Legal Disclaimer

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

To place an order or receive technical assistance

Order/Customer Service: SigmaAldrich.com/order Technical Service: SigmaAldrich.com/techservice Safety-related Information: SigmaAldrich.com/ safetycenter

SigmaAldrich.com

Merck KGaA Frankfurter Strasse 250 64293 Darmstadt, Germany

